
# FastPrep-96™ Pro

## Quick Reference Guide

Revision 2024-11

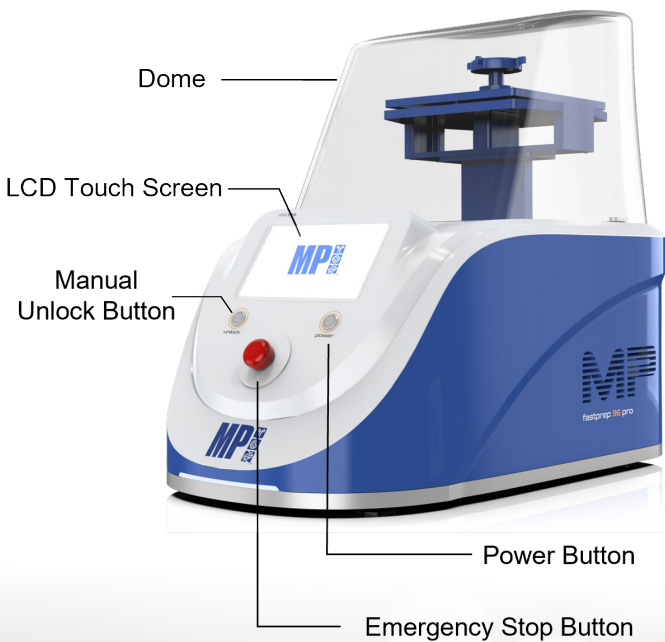### Packing List

Items	SKU	Qty
FastPrep-96™ Pro	116014500	1
Metal QuickFlex™ Adapter	116010575	1
96 x 2 mL tubes		
Blue Clickmaster, Ratchet Nut	116004525	1
Power Cord with Universal Plug	-	1

## Installation & Setup

- Press the unlock button to open the dome manually.

- Remove the foam below the adapter plate.


- Connect the power plug.

- Turn on the power switch located at the back of the machine.

- Power on the machine by pressing the power button located on the front panel.

## Basic Operation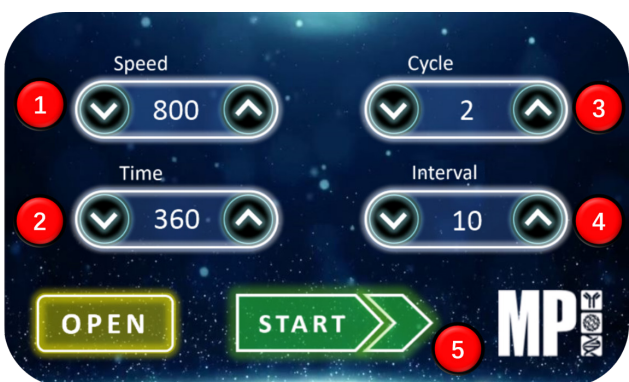

- Press or hold to adjust the speed.
- Press or hold to set the run time in seconds.
- Press to change the number of cycles.
- Press to adjust the duration of the pause between cycles, in seconds.
- Press to start the operation.


REGISTER YOUR INSTRUMENT to activate your warranty

Scan the code or visit [www.mpbio.com/registermyfastprep](http://www.mpbio.com/registermyfastprep)

## | Displays


### Main Screen:

The machine's operational parameters can be configured before starting.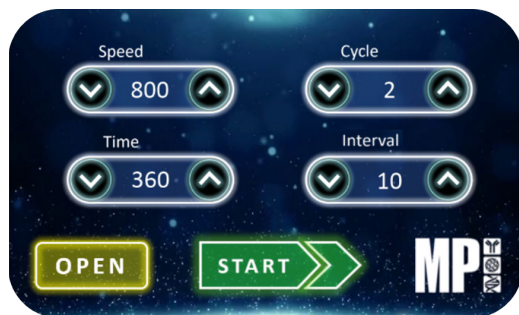


### Running Screen:

The elapsed time is displayed, and the operation can be aborted by pressing the 'STOP' button.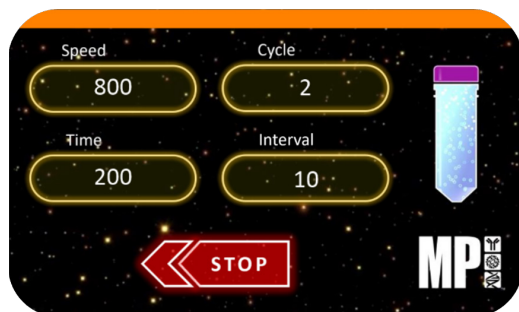

### Cycle Screen:

The interval time is displayed, the operation can be aborted by pressing the 'STOP' button.

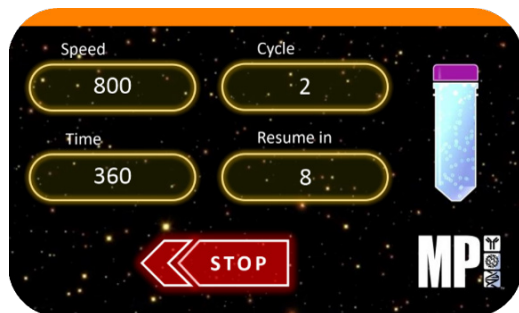## | Safety Features

### Emergency Stop Screen:

This warning screen displays when the Emergency Button is pressed. Twist the knob to reset the machine.

### Dome Opened:

This warning screen displays if the dome is not securely closed before operation.

### For More Information:

Scan this QR code for the instruction manual and video.

## | Accessories List

Metal QuickFlex™ Adapter for 96 x 2 mL tubes for FastPrep-96™ Pro  
(SKU: 116010575)

Metal Dual Plate Holder Assembly for FastPrep-96™ Pro  
(SKU: 119696169)

Blue Clickmaster, Ratchet Nut for FastPrep-96™ Pro  
(SKU: 116004525)

## | Helpful Tips

Wipe the instrument with a damp, soft cloth, and a non-corrosive cleaning agent.

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