



ДИА-М
современная лаборатория

www.dia-m.ru
sales on-line

XPZ PRODUCT MANUALS

TECHNOLOGY LEADING THE INDUSTRY
QUALITY DECIDES THE MARKET





Stablize Fast Safe



CONTENT

01

ABOUT US

02

PRODUCT
FEATURES

03

LABORATORY
GLASSWARE WASHER

04

SPECIAL
CLEANING BASKET

05

BASKET
CATALOGUE



About US

XPZ is a leading manufacturer of laboratory glassware washer, located in Hangzhou city, Zhejiang Province, China. XPZ specializes in research, production and trade the automatic glassware washer which is applied to Bio-pharma, Medical health, Quality inspection environment, Food monitoring, and Petrochemical field.

XPZ is committed to help solve all kinds of cleaning problems. We are the main supplier to Chinese inspection authorities and chemical enterprises, meanwhile, XPZ brand has been spread to many other countries, like India, UK, Russia, Africa, some Europe countries etc., XPZ provides integrated solutions based on customized demand, including product selection, installation and operate training etc.



Passed the
certification
system



We will gather more enterprise advantage to provide innovative products with high quality and excellent service, to keep our long-term friendship.



Modularization




+ Module Frame

Each floor can put two independent module basket racks. One or more floors can be selected for cleaning without changing the module frames.

+ Module

Modules can be freely matched, users can do mix and match according different demands to achieve the cleaning purpose.

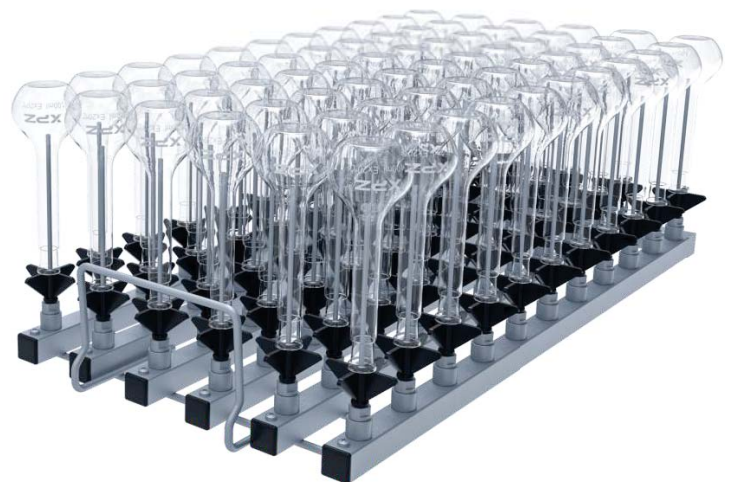

Operation interface

Waterproof stainless steel metal keyboard



OLED color display, spontaneous light, high contrast, no visual angle limit.



+ Basket identification function

Save water, save electricity
Save detergent



+ Password protection

Multi-level password protection



+ Conductivity monitoring

Automatic monitoring water quality



+ Wash start delay function

Time start
Timer start



+ Language control

Support Chinese, English and Russian
three languages



+ Power-off memory function


Can continue wash or cancel wash

Drying System/ Visual viewing window



+ Hot air drying system

Drying loading: Double layer HEPA filter cotton, to further ensure the air purity in the drying process, the instrument with HEPA filter cotton has failure alarm warning, drying time and temperature can be adjusted.

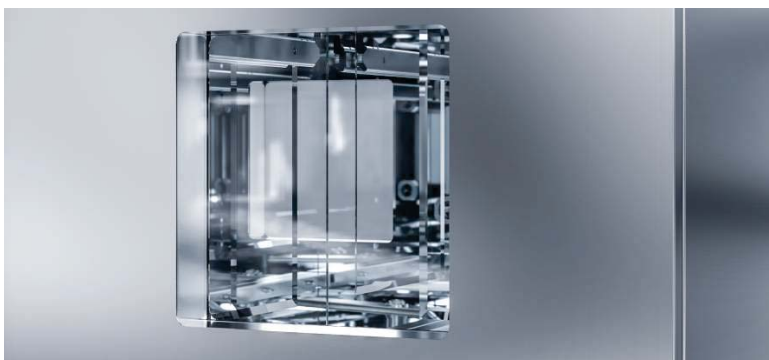


+ Visual viewing window

A large perspective window provides a clear view of the running state.

Double layer toughened glass has good sound insulation and heat insulation effect, with clear and precise texture.

The dual light control makes the entire cleaning process visible.

Automatic induction door technology

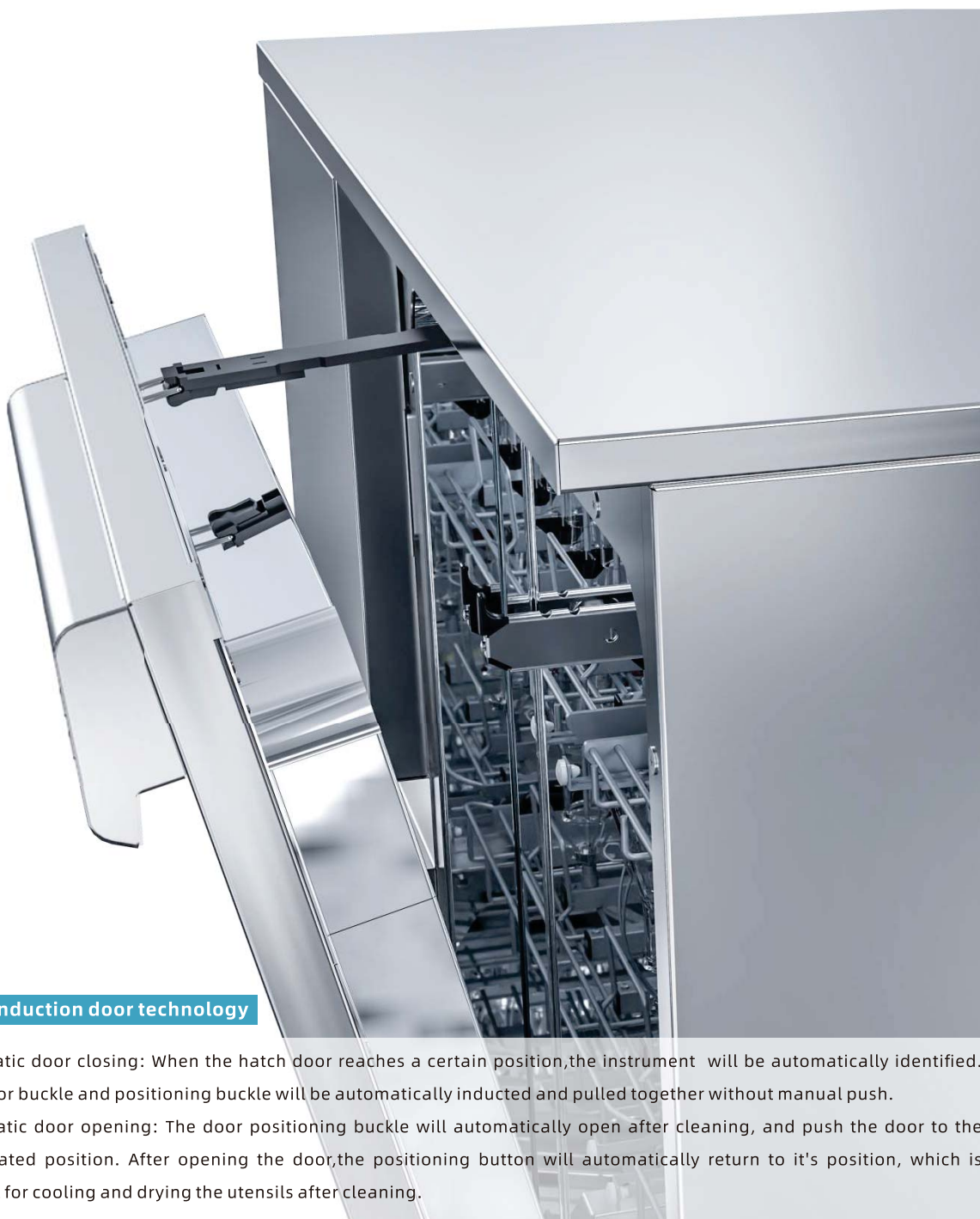

+ ITL induction door technology

Automatic door closing: When the hatch door reaches a certain position, the instrument will be automatically identified. The door buckle and positioning buckle will be automatically inducted and pulled together without manual push.

Automatic door opening: The door positioning buckle will automatically open after cleaning, and push the door to the designated position. After opening the door, the positioning button will automatically return to its position, which is helpful for cooling and drying the utensils after cleaning.

Moment Series

+ Benchtop Style

Moment series are used for cleaning small size utensils.



Moment-F2

Specifications:

Series	Item	Moment-2	Moment-F2
Moment	Dimension	W612*D750*H685mm +Feet10-24mm adjustable	W612*D750*H685mm +Feet10-24mm adjustable
	Chamber volume	126L(Built-in light control)	126L(Built-in light control)
	Number of cleaning layers	1-2layers	1-2layers
	Circulation pump flow rate	320L/min	320L/min
	Heating Power	4kw/9kw	4kw/9kw
	Electricity	220V/380V	220V/380V
	Drying way	Ecodry	Hot air drying
	Instllation method	Benchtop	Benchtop



Feature Description

Series	Item	Moment-2/ Moment-F2
Moment	Operation interface	Microcomputer intelligent control.OLED display, water proof Stainless steel metal buttons with 3 shortcut program keys Support Chinese/Russian/English three languages
	Programs	35 Default,100 user-defined
	Password	Multi-level password protection
	Door	Automatic opening door with viewing window
	Baskets	Modularity basket, one layer can put 2 baskets
	Detection& tracking system	Conductivity detection function(Optional); Printingfunction(Optional); Data storage function
	Wash start delay	Wash start delay function & Delay shutdown function

Glory Series

+ Undercounter Style


Glory series can put under the laboratory table.



Glory-F2

Specifications:

Series	Item	Glory-2	Glory-F2
Glory	Dimension	W612*D750*H830mm +Feet10-24mm adjustable	W612*D750*H830mm +Feet10-24mm adjustable
	Chamber volume	170L(Built-in light control)	170L(Built-in light control)
	Number of cleaning layers	1-3layers	1-3layers
	Circulation pump flow rate	500L/min	500L/min
	Heating Power	4kw/9kw	4kw/9kw
	Electricity	220V/380V	220V/380V
	Drying way	Ecodry	Hot air drying
	Instllation method	Undercounter / Freestanding	Undercounter / Freestanding

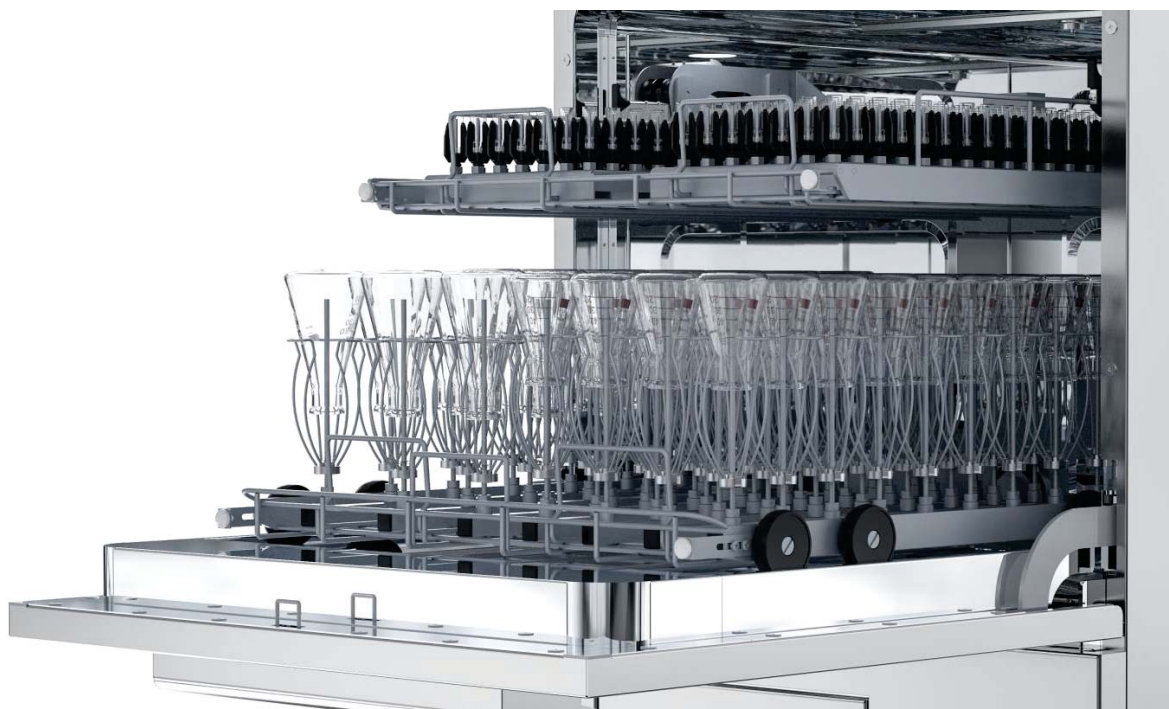

Feature Description

Series	Item	Glory-2/ Glory-F2
Glory	Operation interface	Microcomputer intelligent control.OLED display, water proof Stainless steel metal buttons with 3 shortcut program keys Support Chinese/Russian/English three languages
	Programs	35 Default,100 user-defined
	Password	Multi-level password protection
	Door	Automatic opening door with viewing window
	Baskets	Modularity basket, one layer can put 2 baskets
	Detection& tracking system	Conductivity detection function(Optional); Printingfunction(Optional); Data storage function
	Wash start delay	Wash start delay function & Delay shutdown function
	Basket identification function	Can save water, save electricity, save detergent

Aurora Series

+ Freestanding Style

Aurora series is a freestanding style washer.



Aurora-2/Aurora-F2 PLUS

Specifications:

Series	Item	Aurora-2	Aurora-F2	Aurora-F2 PLUS
Aurora	Dimension	W612*D750*H990mm +Feet10-24mm adjustable	W930*D750*H990mm +Feet10-24mm adjustable	W930*D750*H1110mm +Feet10-24mm adjustable
	Chamber volume	202L(Built-in light control)	202L(Built-in light control)	239L(Built-in light control)
	Number of cleaning layers	1-3layers	1-3layers	1-3layers
	Circulation pump flow rate	0-600L/min	0-600L/min	0-600L/min
	Heating Power	4kw/9kw	4kw/9kw	4kw/9kw
	Electricity	220V/380V	220V/380V	220V/380V
	Drying way	Ecody	Hot air drying	Hot air drying
	Instllation method	Freestanding	Freestanding	Freestanding



Aurora-F2

Feature Description

Series	Item	Aurora-2/Aurora-F2/Aurora-F2 PLUS
Aurora	Operation interface	Microcomputer intelligent control.OLED display, water proof Stainless steel metal buttons with 3 shortcut program keys Support Chinese/Russian/English three languages
	Programs	35 Default,100 user-defined
	Password	Multi-level password protection
	Door	Automatic opening door with viewing window
	Baskets	Modularity basket, one layer can put 2 baskets
	Detection& tracking system	Conductivity detection function(Optional); Printingfunction(Optional); Data storage function
	Wash start delay	Wash start delay function & Delay shutdown function
	Basket identification function	Can save water, save electricity, save detergent

Flash Series

+ Freestanding Style

Flash series is a freestanding style washer.



Flash-F2

Specifications:

Series	Item	Flash-F2	Flash-F2 PLUS
Flash	Dimension	W930*D750*H1326mm +Feet 82-90mm adjustable	W930*D750*H1476mm +Feet 82-90mm adjustable
	Chamber volume	308L(Built-in light control)	345L(Built-in light control)
	Number of cleaning layers	1-4layers	1-4layers
	Circulation pump flow rate	0-1000L/min	0-1000L/min
	Heating Power	4kw /9kw / 18kw	4kw /9kw / 18kw
	Electricity	220V/380V	220V/380V
	Drying way	Hot air drying	Hot air drying
	Instllation method	Freestanding	Freestanding



Flash-F2 PLUS

Feature Description

Series	Item	Flash-F2/Flash-F2 PLUS
Flash	Operation interface	Microcomputer intelligent control.OLED display, water proof Stainless steel metal buttons with 3 shortcut program keys Support Chinese/Russian/English three languages
	Programs	35 Default,100 user-defined
	Password	Multi-level password protection
	Door	Automatic opening door with viewing window
	Baskets	Modularity basket, one layer can put 2 baskets
	Detection& tracking system	Conductivity detection function(Optional); Printingfunction(Optional); Data storage function
	Wash start delay	Wash start delay function & Delay shutdown function
	Basket identification function	Can save water, save electricity, save detergent

Module Frame



**Upper or Middle level basket frame
FA-Z01**

- With two connectors
- Used to connect 2 injection baskets
- Auto self-sealing docking valve
- H139, W529, D570mm




**Upper or Middle level basket frame
FA-Z03**

- To load non-injection baskets
- Built-in spray arm
- H183, W530, D569mm



**Upper or Middle level basket frame
FA-Z05**

- To load non-injection baskets
- Built-in spray arm
- H139, W529, D574mm

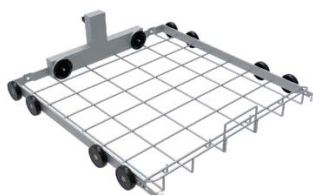

**Lower level basket frame
FA-Z02**

- With two connectors
- Used to connect 2 injection baskets
- Auto self-sealing docking valve
- H148, W531, D577mm




**Lower level basket frame
FA-Z04**

- To load non-injection baskets
- H83, W531, D560mm



**Lower level basket frame
FA-Z06**

- Double levels
- To load non-injection baskets
- Built-in spray arm
- H230, W528, D584mm

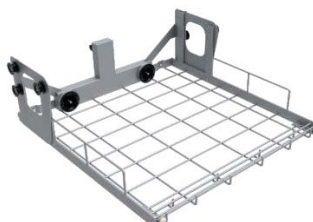

**Lower level basket frame
FA-Z07**

- With two connectors
- Used to connect 2 injection baskets
- Auto self-sealing docking valve
- H182, W529, D564mm



**Full frame injection module
FA-Z11**

- 38 pipettes can be loaded in three layers
- 1st Row:10pcs 10-100ml pipettes,H550mm
- 2nd Row:14pcs 10-25ml pipettes,H500mm
- 3rd Row:14pcs 1-10ml pipettes,H440mm
- H373, W531, D582mm

Cleaning Module



Injection Module 3 injections FA-K03

- For 500-2000ml Erlenmeyer flasks, Reagent bottle, Blue cap bottle, 500-1000ml Volumetric flask and so on.
- Injector nozzle: $\phi 6 \times H230\text{mm}$
- H258, W194, D482mm



Injection Module 9 injections FA-L09

- For 500-2000ml Erlenmeyer flasks, Reagent bottle, Blue cap bottle, 500-1000ml Volumetric flask and so on.
- Injector nozzle: $\phi 6 \times H230\text{mm}$
- H263, W499, D505mm



Injection Module 10 injections FA-K10

- For 100-500ml Erlenmeyer flasks, Reagent bottle, Blue cap bottle, Volumetric flask and so on. 10-100ml cylinder.
- Injector nozzle: $\phi 6 \times H230\text{mm}$
- H249, W140, D482mm



Injection Module 10 injections FA-K11

- For 100-500ml Erlenmeyer flasks, Reagent bottle, Blue cap bottle, Volumetric flask and so on. 10-100ml cylinder.
- Injector nozzle: $\phi 6 \times H230\text{mm}$
- H249, W140, D482mm



Injection Module 10 injections FA-K11/1

- For 250-500ml Erlenmeyer flasks, Reagent bottle, Blue cap bottle, Volumetric flask and so on.
- Injector nozzle: $\phi 6 \times H130\text{mm}$
- H179, W240, D505mm
- Holder's height can not adjustable



Injection Module 12 injections FA-K12/1

- Suitable for $\phi 3.5\text{mm}$ Viscometer
- Injector nozzle: $\phi 2.5 \times H110\text{mm}$
 $\phi 6 \times H230\text{mm}$
- H249, W172, D482mm



Injection Module 14 injections FA-K14

- For 50-250ml Erlenmeyer flasks, Volumetric flask, 10-50ml cylinder and so on.
- Injector nozzle: $\phi 4 \times H160\text{mm}$
- H179, W163, D482mm



Injection Module 15 injections FA-K15

- For 50-250ml Erlenmeyer flasks, Reagent bottle, Blue cap bottle, Volumetric flask and so on.
- Injector nozzle: $\phi 4 \times H160\text{mm}$
- H179, W175, D482mm



Injection Module 18 injections FA-K18

- For 100-250ml Erlenmeyer flasks, Reagent bottle, Blue cap bottle, 10-50ml cylinder and so on.
- Injector nozzle: $\phi 4 \times H140\text{mm}$
- H167, W232, D482mm

Cleaning Module




Injection Module 21 injections FA-K20

- For 50-150ml Erlenmeyer flasks, Reagent bottle, Volumetric flask, 10-50ml cylinder and so on.
- Injector nozzle: $\phi 4 \times H160\text{mm}$
- H179, W192, D482mm



Injection Module 21 injections FA-K21

- For 50-150ml Erlenmeyer flasks, Reagent bottle, Blue cap bottle, Volumetric flask and so on.
- Injector nozzle: $\phi 4 \times H160\text{mm}$
- H179, W192, D482mm

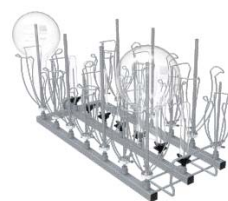

Injection Module 21 injections FA-K22

- For 50-150ml Erlenmeyer flasks, Reagent bottle, Blue cap bottle, Volumetric flask and so on
- Injector nozzle: $\phi 2.5 \times H110\text{mm}$
 $\phi 4 \times H160\text{mm}$ $\phi 6 \times H230\text{mm}$
- H249, W192, D482mm




Injection Module 36 injections FA-K30

- For 25-100ml Erlenmeyer flasks, Reagent bottle, Volumetric flask, 5-10ml test tube and so on.
- Injector nozzle: $\phi 2.5 \times H110\text{mm}$
- H129, W200, D482mm



Injection Module 36 injections FA-K31

- For 25-100ml Erlenmeyer flasks, Reagent bottle, Volumetric flask and so on
- Injector nozzle: $\phi 4 \times H160\text{mm}$
- H179, W200, D482mm

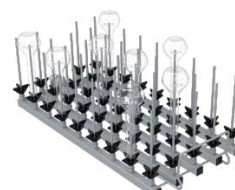

Injection Module 44 injections FA-K44

- For 25-50ml Volumetric flask, test tube, vials and so on
- Injector nozzle: $\phi 4 \times H140\text{mm}$
- H159, W215, D482mm



Injection Module 66 injections FA-K66

- For 25ml Volumetric flask, test tube, vials and so on
- Injector nozzle: $\phi 2.5 \times H110\text{mm}$
- H129, W215, D482mm



Injection Module 119 injections FA-K90

- For 1.5ml vials specially
- Injector nozzle: $\phi 2.5 \times H53\text{mm}$
- H76, W190, D488mm



Injection Module 119 injections FA-K91

- For pipettes specially
- H191, W190, D488mm

Cleaning Module / Basket / Cover



Injection Module 65 injections

FA-M65

- Can load 28pcs pipettes, 28pcs vials, 9pcs Erlenmeyer flasks, volumetric flask, cylinder and so on.
- Injector nozzle: $\phi 2.5 \times H110\text{mm}$; $\phi 4 \times H160\text{mm}$; $\phi 6 \times H230\text{mm}$
- H257, W190, D488mm



Injection Module 112 injections

FA-M112

- Can load 56pcs pipettes, 56pcs test tubes and so on.
- Injector nozzle: $\phi 2.5 \times H110\text{mm}$
- H184, W190, D488mm



Full Basket 56

T-101

- Suitable for clean petri dishes
- 56 seats
- H145, W485, D504mm



FullBasket 56

T-102

- Only used together with T-101
- Suitable for clean petri dishes
- 56 seats
- H145, W485, D504mm



Full Basket

T-103

- For funnel, beakers, jar, etc
- H120, W446, D470mm

Cover G-204

- The cover for T-103
- H28, W223, D423mm



Cover net

G-201/G-202


- Basket covering for bottles to prevent the glassware rushed out.
- G-201 is suitable for FA-K90
- G-202 is suitable for T-101/T-102
- H10, W215, D455mm



Shelf

T-203

- 21 spring clips are used to hold the utensils
- The distance between the clips is about 60mm
- H155, W230, D460mm

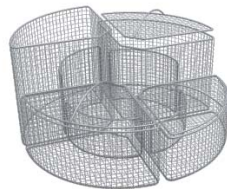

1/4 Basket

T-401/402/403/404

- H100, W218, D218mm, The maximum is 12*75mm
- H127, W218, D218mm, The maximum is 12*105mm
- H187, W218, D218mm, The maximum is 12*165mm
- H230, W218, D218mm, The maximum is 12*200mm



Half Basket

T-204/2


- For funnel, beakers, jar, etc
- H119, W224, D440mm

Cover G-401

- The cover for T-204/2
- H21, W210, D210mm

Laboratory Glassware Washer

● Means support this model
○ Means not support this model

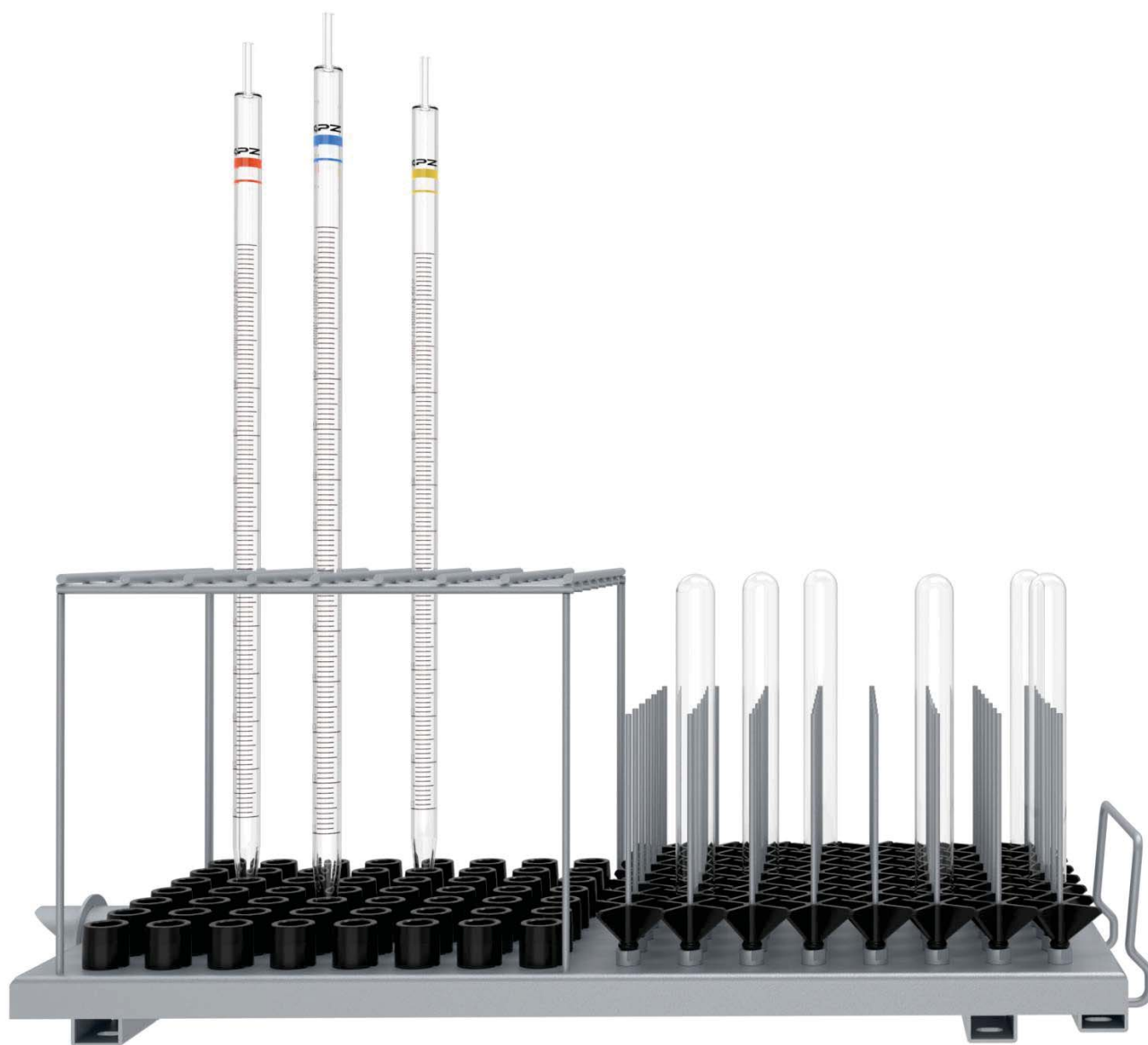

Only do professional service for you

000 «Диаэм»

Москва

ул. Магаданская, д. 7, к. 3 ■ тел./факс: (495) 745-0508 ■ sales@dia-m.ru

www.dia-m.ru

С.-Петербург

+7 (812) 372-6040
spb@dia-m.ru

Новосибирск

+7 (383) 328-0048
nsk@dia-m.ru

Воронеж

+7 (473) 232-4412
vrn@dia-m.ru

Йошкар-Ола

+7 (927) 880-3676
nba@dia-m.ru

Красноярск

+7 (923) 303-0152
krsk@dia-m.ru

Казань

+7 (843) 210-2080
kazan@dia-m.ru

Ростов-на-Дону

+7 (863) 303-5500
rnd@dia-m.ru

Екатеринбург

+7 (912) 658-7606
ekb@dia-m.ru

Кемерово

+7 (923) 158-6753
kemerovo@dia-m.ru

Самара

+7 (915) 482-4897
samara@dia-m.ru

