

Instructions for use

Eppendorf Heat Sealing Film & Foil

English (EN)

Eppendorf Heat Sealing Film & Foil

1 Product description

Eppendorf Heat Sealing Film and Foil are designed for sealing PCR plates as well as micro test plates and deepwell plates made of polypropylene. They provide optimum protection against evaporation when used in conjunction with a suitable heat sealer. They are suitable for long term storage.

The transparent Heat Sealing Film allows sample monitoring and is suitable for fluorescence measurements. The Heat Sealing Foil can be easily pierced by pipette tips and can be removed again from the plate.

The products are designed for single use in routine, research and training labs in the bioscience, industrial and chemical fields. They are intended for use by trained specialists only.

Adhesive edges with white cover strips on the underside.



2 Safety





WARNING! Damage to health from toxic, radioactive or aggressive chemicals as well as infectious liquids and pathogenic germs.

- ▶ Observe the national regulations for handling these substances, the biological security level of your laboratory, the material safety data sheets and the manufacturer's application notes.
 - ▶ Wear personal protective equipment.
 - ▶ For full instructions regarding the handling of germs or biological material of risk group II or higher, please refer to the "Laboratory Biosafety Manual" (Source: World Health Organization, current edition of the Laboratory Biosafety Manual).
-
- ▶ Make sure to read the operating manuals of the respective heat sealers and plates prior to use. Note in particular the high temperatures required for sealing films and foils.
 - ▶ Reagents, disposables contaminated with reagents and materials used for cleaning and disinfecting should be disposed of in accordance with laboratory regulations.
 - ▶ Use only undamaged articles.
 - ▶ Use articles from undamaged packages only.

3 Application

1. Use gloves.
2. Peel off the cover strips.
3. Position the film or the foil with the adhesive edges on the plate.
4. Seal the plate using the heat sealer. The optimum time depends on the plate type. Start with 2 to 3 s.

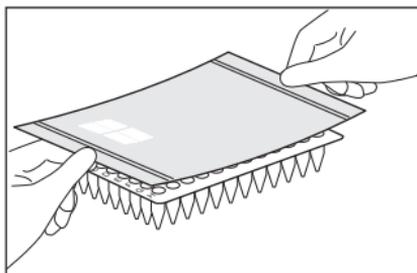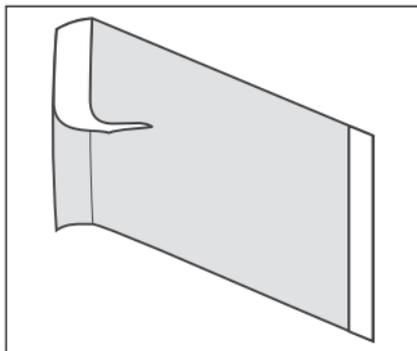

4 Technical data

Eppendorf PCR clean	Certified free of RNase, DNase, human DNA and PCR inhibitors*.
Storage	Protect from direct sunlight and UV light. Store dry at room temperature.
Operational temperature	-80°C to +110°C

* Lot-specific certificates are available at www.eppendorf.com or certificates@eppendorf.de.

5 Ordering information

Order no. (International)	Order no. (North America)	Description
0030 127.838	0030127838	Heat Sealing Film 100 pieces
0030 127.854	0030127854	Heat Sealing Foil 100 pieces
5391 000.001	5391000001	Heat Sealer 110 V – 230 V / 50 Hz– 60 Hz S100
5392 000.005	5392000005	S200

000 «Диаэм»

Москва
ул. Магаданская, д. 7, к. 3 ■ тел./факс: (495) 745-0508 ■ sales@dia-m.ru

www.dia-m.ru

С.-Петербург
+7 (812) 372-6040
spb@dia-m.ru

Новосибирск
+7(383) 328-0048
nsk@dia-m.ru

Воронеж
+7 (473) 232-4412
vrn@dia-m.ru

Йошкар-Ола
+7 (927) 880-3676
nba@dia-m.ru

Красноярск
+7(923) 303-0152
krsk@dia-m.ru

Казань
+7(843) 210-2080
kazan@dia-m.ru

Ростов-на-Дону
+7 (863) 303-5500
rnd@dia-m.ru

Екатеринбург
+7 (912) 658-7606
ekb@dia-m.ru

Кемерово
+7 (923) 158-6753
kemerovo@dia-m.ru

Армения
+7 (094) 01-0173
armenia@dia-m.ru

